

## PAL RTC Grabber DW-96

### Automated Transport of 96 well plates ...


- Transport of 96 well plates
- Automatic solvent transfer to 96 well plates
- Shaker up to 5'000 rpm
- Optional Heater

### You need a special solution? Challenge us ....

- Send us your request and we will make it possible




Add solvent to 96 well plate

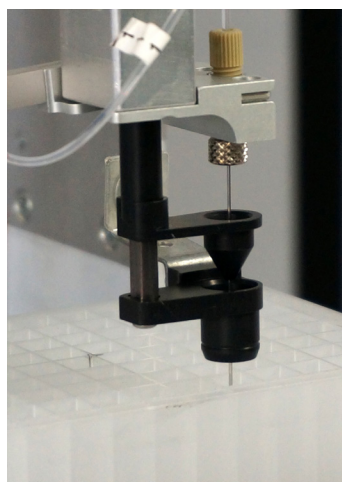

Pick-up Grabber DW-96



Place 96 well plate on shaker or centrifuge

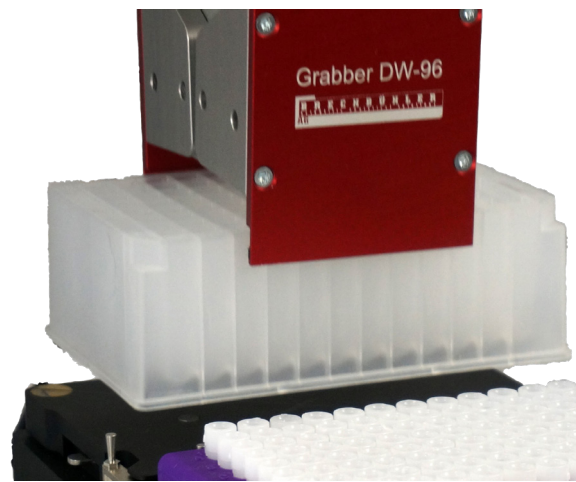


### Automated Transport of anything you can think of ...

- Transport of any item fitting to the 96 well plates format
- Customized workflows available on request
- Systems delivered with standard tasks
- Options available:  
96 well plates Centrifugation up to 2'000 g  
Heated shaker  
Sonication
- All PAL options available such as:  
Bar code reader  
Injection into LC (LC/MS) system  
And much more


## PAL RTC Grabber DW-96

### DW-96 starter kit ...


- Grabber D885-SL
- Trayholder legs 287 mm (set of 2)    Pos. 1
- Grabber DW-96    Pos. 2
- Centering Plate DW-96    Pos. 3
- Optional Shaker    Pos. 4

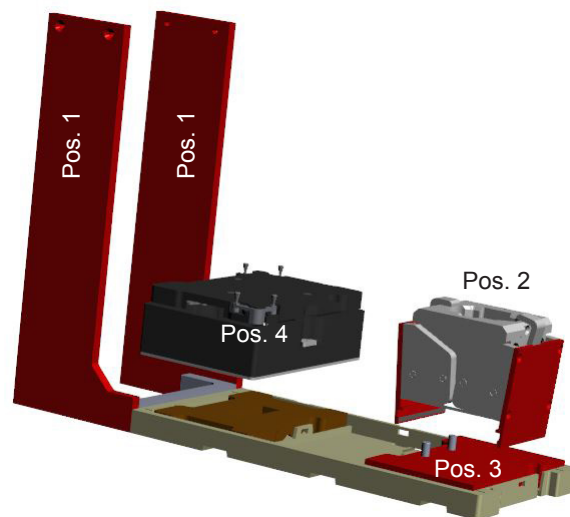

Centering Plate DW-96

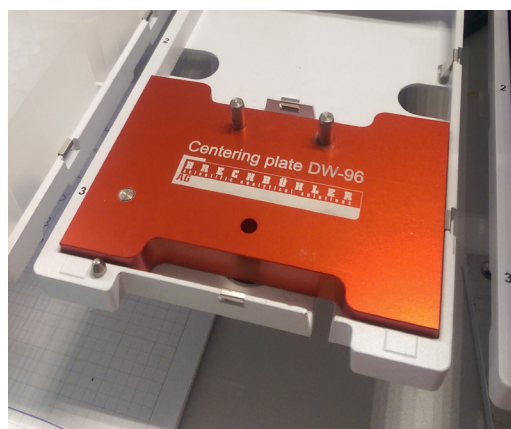

Grabber DW-96 on Centering Plate

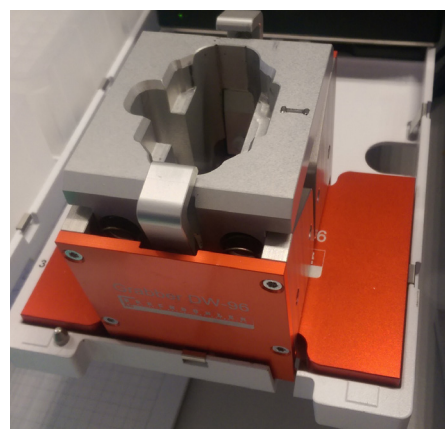

### Method / Tasks provided...

- **Get DW96 grabber**  
Get the D885-SL grabber and get the DW96 grabber
- **Drop DW96 Grabber**  
Return the DW96 grabber to its position
- **Transport plate**  
Transport a plate from a defined position to another one
- **Home shaker**  
Home shaker after a movement
- **Unlock shaker**  
Open the shaker to place a plate
- **Lock shaker**  
Lock a plate on the shaker
- **Start shaker at set speed for set time**  
Start the shaker at the set speed and time
- **Shaker operation requires Brechbuehler control software**  
(included with shaker)